



## Configuring ibidi $\mu$ -Slides

Ibidi provides  $\mu$ -slides with 2-18 wells, which all fit into the sample carrier insert "2x Slide – Long" (for 2 standard slides, 76 x 26 mm). The following procedure shows how to configure ZEN to use a template with a different number of wells.

1. In the "Sample" tab, go to "Prescan Options". Check "Sample Carrier Detection" and click on "Configure...".

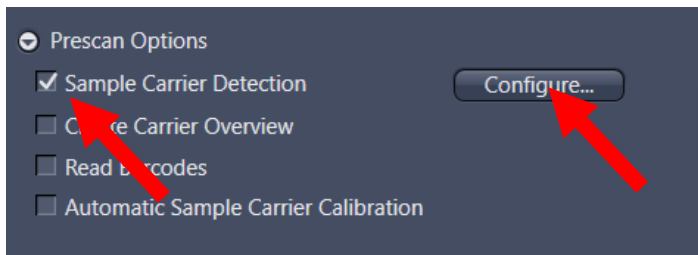

2. In the "Detection Result" list of the dialog, select "Insert – 2x Slide – Long", then click "Select..."

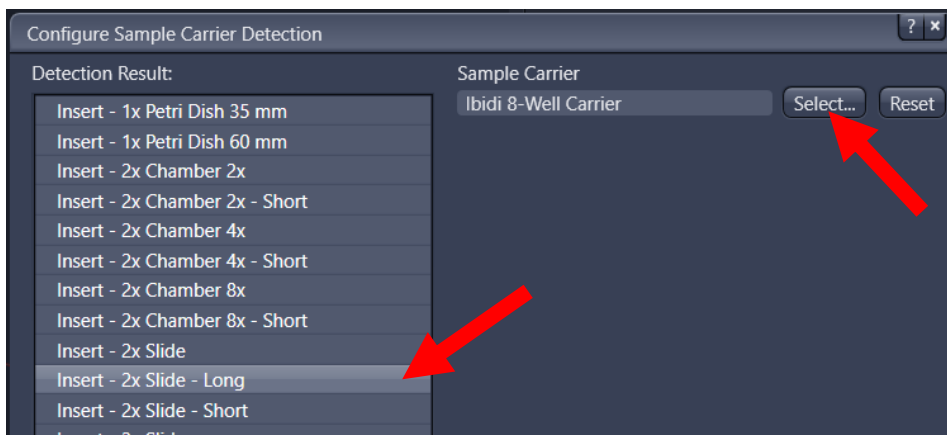

3. In the "Select Template" dialog, select the desired template, then click "OK"

